



 *Formaway*
FORMALIN NEUTRALIZER PRODUCTS

Formaway Formalin Neutralizing Absorbent Pads

Mopec's Formaway Formalin Neutralizing Absorbent Pad (F-NAP) was designed to reduce formaldehyde vapors while dissecting tissue, transferring tissue from containers of formalin, and transporting or storing tissue in containers of formalin, by absorbing and neutralizing formalin. Available in miniature, small, medium, and large. Sold by the pack or by the case.



SKU	Description	Dimensions	Quantity	Absorbancy
BE092	Formalin Neutralizing Pads, Mini	3" x 4"	100/pack	12-13ml/pad
BE093	Formalin Neutralizing Pads, Small	8" x 11"	25/pack	100ml/pad
BE094	Formalin Neutralizing Pads, Medium	11" x 15"	25/pack	200ml/pad
BE095	Formalin Neutralizing Pads, Large	15" x 20"	25/pack	300ml/pad

Formaway Formalin Neutralizing Powder

Mopec's Formaway Formalin Neutralizing Powder safely and easily neutralizes spent formalin waste to a non-hazardous and non-toxic solution in only 20 minutes. Contents are emptied directly into 1 gallon of spent 10% Formalin or 4% Glutaldehyde, reducing risk exposure to lab personnel. This unique formula also eliminates drain-clogging polymers. Available in a case of 12 or 32 bottles.



SKU	Description	Size	Quantity
BG550-12	Formalin Neutralizing Powder	16 oz. bottle	12/case
BG550-32	Formalin Neutralizing Powder	16 oz. bottle	32/case
BG551	Waste Collection Container	1 gallon	each
BG552	Waste Collection Container	2.5 gallons	each
BG553	Funnel with Tissue Screen	-	each
BG554	Formalin Test Strips	-	100/pack

Formaway Green Liquid Formalin Neutralizer

Mopec's Formaway Green Liquid Formalin Neutralizer quickly and effectively neutralizes spent formalin for sewer disposal. Simply add 32 ounces of Formaway Green to each gallon of 10% neutral buffered formalin waste and allow to stand for at least 4 hours before disposing. Available in 1 gallon, 2.5 gallon, and 5 gallon cubes.

SKU	Description	Capacity
BG560-1	Formalin Green Liquid Neutralizer	1 gallon cube
BE560-2	Formalin Green Liquid Neutralizer	2.5 gallon cube
BE560-5	Formalin Green Liquid Neutralizer	5 gallon cube



SCAN THE QR CODE OR VISIT [MOPEC.COM/FORMAWAY](https://www.mopec.com/formaway) FOR MORE INFORMATION