



| REV | DESCRIPTION | DATE | ECN # | APPR |
|-----|------------------|-----------|-------|------|
| A | INITIAL RELEASE | 00-XXX-19 | XXXX | MM |
| B | UPDATE NEW MODEL | XXXX | XXXX | JJ |
| | | | | |
| | | | | |
| | | | | |

| Item # | Qty. | File Name | Title | Material | Rev. |
|--------|------|-----------|--|--------------------------|------|
| 1 | 1 | 017-1 | BAR CODE ASY PURCHASED PART | | A |
| 2 | 1 | 017-2 | BAR CODE ASY PURCHASED PART | | A |
| 3 | 1 | 017-3 | BAR CODE ASY PURCHASED PART | | A |
| 4 | 1 | 017-4 | BAR CODE ASY PURCHASED PART | | A |
| 5 | 2 | PB0063 | WELD STUD - SS, 8-32 x 1/2"L | HYBROCO # CD50832-0500-1 | A |
| 6 | 2 | PB0068 | NUT - SS, 8-32, ACORN NUT | McMASTER-CARR # 91855A44 | A |
| 7 | 2 | PB0156 | NUT - ZINC, 8-32, STANDARD W/ TOOTH WASHER | | A |
| 8 | 1 | S6800-A | PEGBOARD MEDIUM MOUNTING TAB | 18 GA. TYPE T304 SS | A |

THIS DOCUMENT CONTAINS PROPRIETARY INFORMATION OF MOPEC, AND MAY NOT BE COPIED IN WHOLE OR PART, AND NO USE SHALL BE MADE OF IT OTHER THAN FOR WHICH IT HAS BEEN SUBMITTED WITHOUT THE WRITTEN PERMISSION OF A MOPEC. MANAGER

MOPEC 2024

UNLESS OTHERWISE SPECIFIED:
ALL DIMENSIONS IN INCHES

| | | |
|------------|--------|-------------|
| TOLERANCES | 0.0 | ± 0.1 in |
| | 0.00 | ± 0.05 in |
| | 0.000 | ± 0.010 in |
| | 0.0000 | ± 0.0010 in |
| ANGLES | | ± 0.5° |

BREAK ALL CORNERS 0.063 X 45°
ALL FILLET RADII R0.063 /0.125in
INSPECT OVAL DIMENSIONS

DRAWN TO 3RD ANGLE PROJECTION

MOPEC GROUP

MOPEC MORTTECH

ADVANCED TECHNOLOGY MANUFACTURING

DRAWN BY: B. LAY
DATE: 2/20/2023

CHECKED BY: ---
DATE: 2/20/2023

APPROVED BY: M. McCLAIN
DATE: 2/20/2023

PART NUMBER: AX017-B

DESCRIPTION: BARCODE HOLDER

SIZE: A2
FILE NAME: AX017-B.dft
REV: B

SCALE: 1:1

DO NOT SCALE IF IN DOUBT ASK

SHEET 2 OF 2